

KEYWORDS:

Herbicides

Triazines


Notes: BPX35 provides quick analysis of all 3 Triazine compounds

HERBICIDES

Part No.: 054711
Phase: BPX35, 0.25µm film
Column: 25m x 0.22mm ID
Initial Temp: 80°C
Rate : 10°C/min
Final Temp : 300°C 5min
Detector: FID, 380°C
Injection Mode: Split (20:1)
Carrier Gas: He, 100 kpa

Components

- | | |
|-------------------------|--------------------------|
| 1. Eptam [®] | 12. Atrazine |
| 2. Sutan [®] | 13. Simazine |
| 3. Vernam [®] | 14. Terbacil |
| 4. Tillam [®] | 15. Sencor [®] |
| 5. Ordram [®] | 16. Dual [®] |
| 6. Treflan [®] | 17. Paarlan [®] |
| 7. Balan [®] | 18. Prowl [®] |
| 8. Ro-Neet [®] | 19. Bromacil |
| 9. Propachlor | 20. Oxadiazon |
| 10. Tolban [®] | 21. GOAL [®] |
| 11. Propazine | 22. Hexazinone |

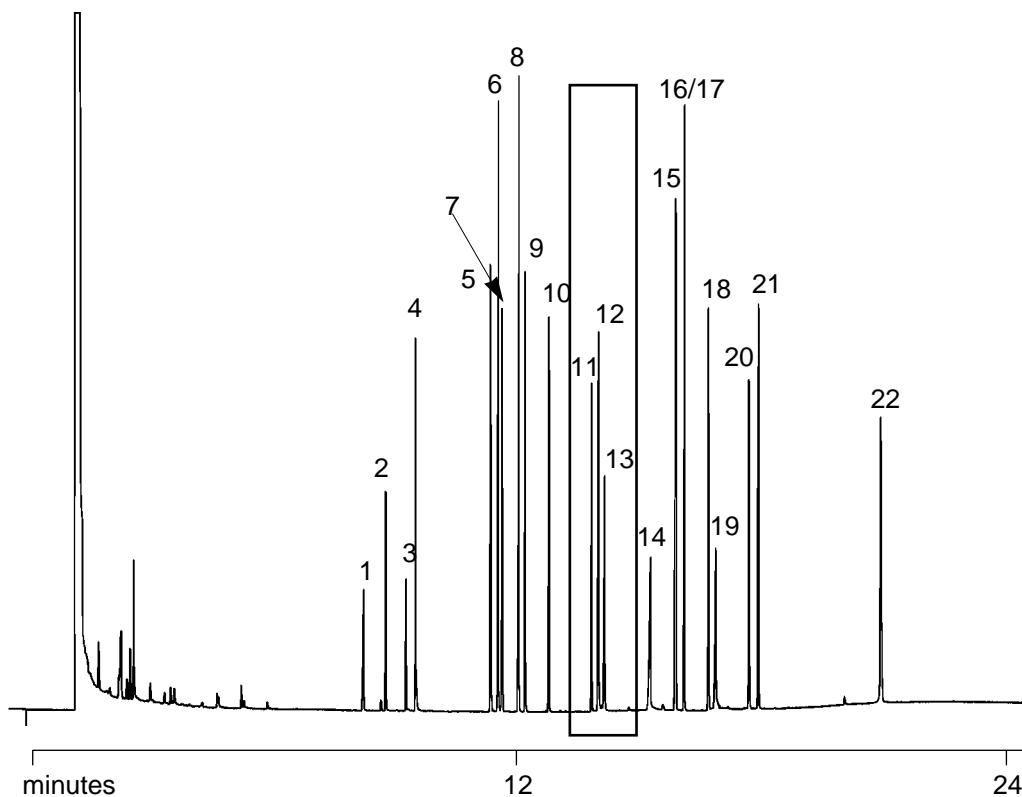

visit www.sge.com or contact our technical Customer Support Team on: techsupport@sge.com

SGE International Pty. Ltd.
email: support@sge.com

SGE Incorporated (USA)
email: usa@sge.com

SGE (France) Sarl
email: france@sge.com

SGE Deutschland GmbH
email: germany@sge.com

SGE Japan Inc.
email: sgejpn@itjit.ne.jp

SGE Australia Pty. Ltd.
email: sales@sge.com

SGE Europe Ltd. (UK)
email: uk@sge.com

SGE (Italia) Srl.
email: sge.italia@tin.it

SGE China Service Centre
email: sge@bj.col.com.cn

SGE India
email: sgeindia@bom4.vsnl.net.in